

What you can't see can't hurt you / Millipore Corporation.

Contributors

Millipore Corporation.

Publication/Creation

Bedford, MA : Millipore, 1995.

Persistent URL


<https://wellcomecollection.org/works/rduy4hxc>

License and attribution

Conditions of use: it is possible this item is protected by copyright and/or related rights. You are free to use this item in any way that is permitted by the copyright and related rights legislation that applies to your use. For other uses you need to obtain permission from the rights-holder(s).



Wellcome Collection
183 Euston Road
London NW1 2BE UK
T +44 (0)20 7611 8722
E library@wellcomecollection.org
<https://wellcomecollection.org>

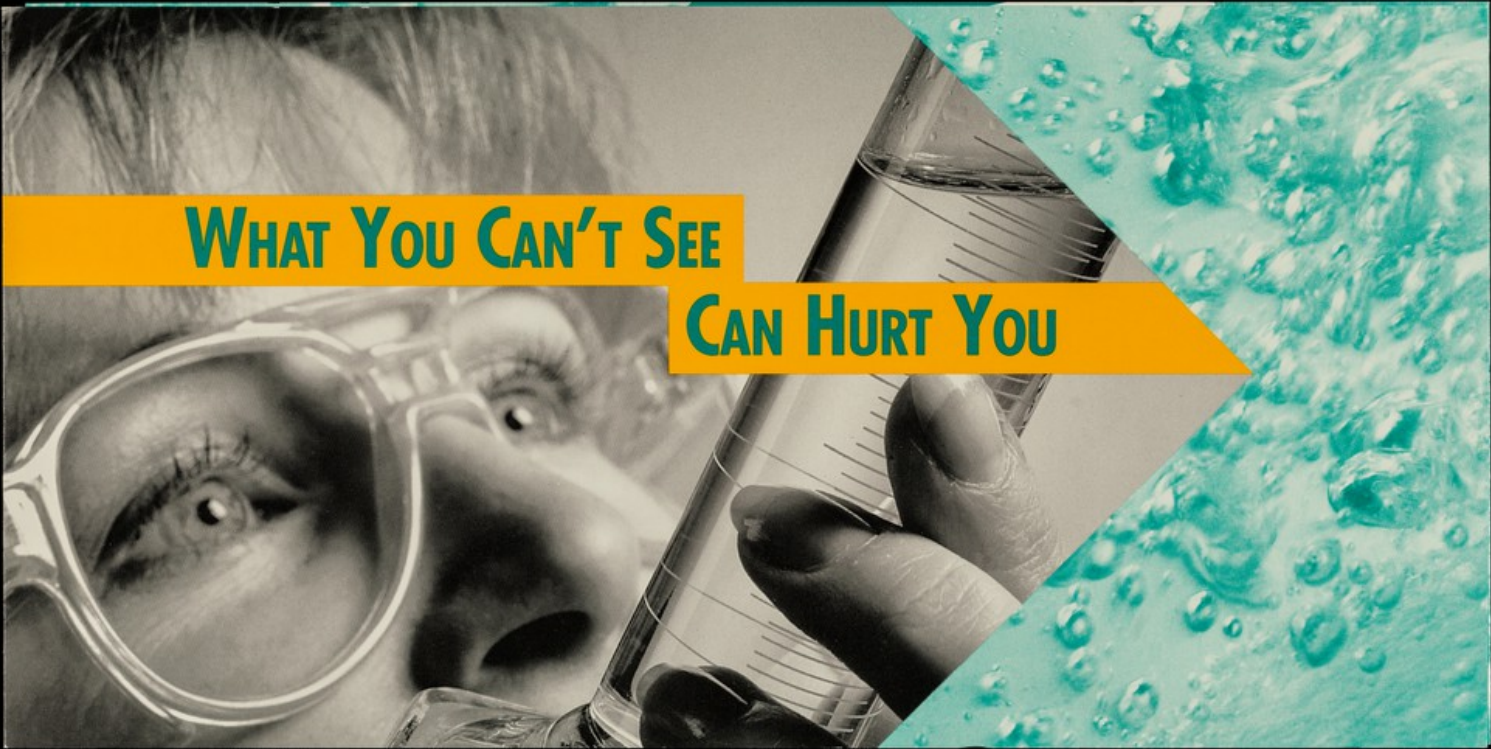


WHAT YOU CAN'T SEE

CAN HURT YOU

Do You Know What's Really In Your Ultrapure Water?

When it comes to ultrapure water, what you can't see can often hurt your application. Your system's resistivity meter may indicate that the water produced is at the theoretically "pure" limit of 18.2 megohm-cm, but don't be fooled! Resistivity is only a measure of ionic purity. In spite of this maximum resistivity reading, your water can still contain high concentrations of organic contaminants that go undetected by your system's resistivity monitor.

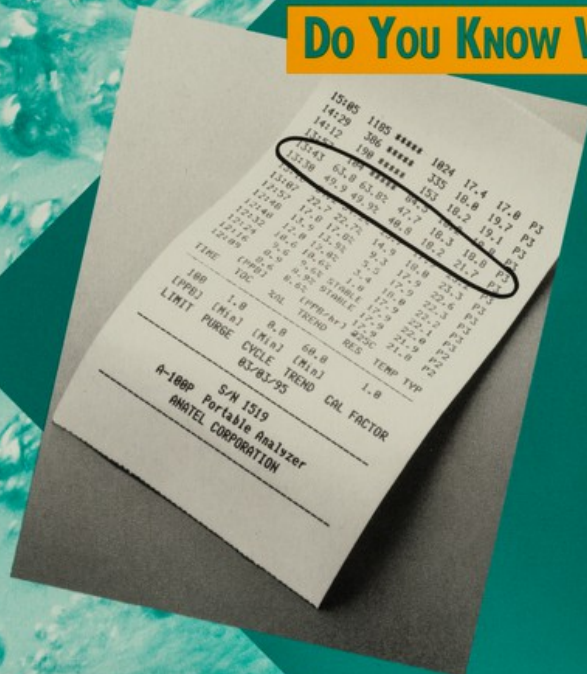

**Water purity you can see
with the new Millipore
A-10 TOC Monitor**

Now you can be sure of organic water purity with the new Millipore Anatel A-10 TOC Monitor. It can be simply connected to your water purification system to provide quantitative measurement of organic purity.

The A-10 monitor's single mode key allows reporting of all data, and operating conditions simply and conveniently. And the LCD displays the values and parameters in a large, easy-to-see format.



Millipore now offers the A-10 TOC monitor developed with Anatel Corporation, the world leader in TOC measurement for high purity water production in the microelectronics industry.

**Now see the
complete picture of
water purity**

When used with the **Milli-Q[®] 185 Plus Ultrapure Water System**, the A-10 monitor assures you of ultra-low organics water purity you can see for your critical organic-sensitive analyses.

The Milli-Q[®] 185 Plus system represents a new advance in water purity, utilizing UV photo-oxidation to provide water with extremely low TOC levels for applications such as environmental monitoring, HPLC, IC, GC, and toxicology studies.



**Get results
you can actually see**


Don't stay in the dark. Get results you can actually see. To learn more about the Milli-Q[®] 185 Plus system, including our product brochure and technical brief, as well as the new Anatel A-10 monitor, call our technical specialists at **01923 816375**, or fax **01923 816128**.

Measuring resistivity is not enough

It's a fact that most ultrapure water purification systems exhaust their capacity to remove dissolved organic contaminants before they lose their capacity to remove ions. The resulting "organics breakthrough" leads to high and variable organic levels in the product water.

With today's demanding instrumentation, even low concentrations of dissolved organic compounds in ultrapure water can interfere with analyses and ruin your results. Not only do you require a system which provides ultra-low organics water, but you also need a way to measure organic purity.

Organic Breakthrough

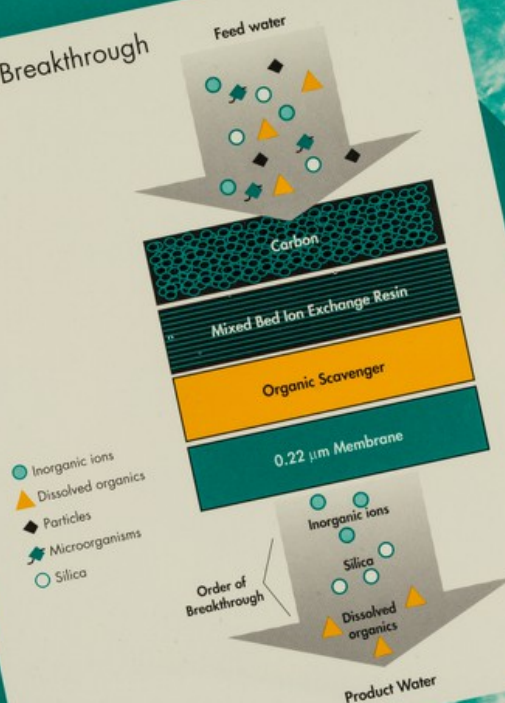

MILLIPORE

$(H_2O)^\infty$

Millipore is a registered trademark of Millipore Corporation or an affiliated company.

Milli-Q is a trademark of Millipore Corporation or an affiliated company.

Anatel is a trademark of Anatel Corporation.

Lit. No. DM022 Printed in U.S.A. 5/95 95-200 UK

Copyright © 1995 Millipore Corporation or an affiliated company, Bedford, MA

Austria, Central Europe, C.I.S., Africa, Middle East and Gulf: 43 (1) 877 8926

Belgium & Luxembourg: (02) 726 88 40 **Denmark:** 46 59 00 23

Finland: (90) 804 51 10 **France:** (1) 30 12 70 00

Germany: (06196) 494-0 **Italy:** (02) 25078 1

The Netherlands: (01608) 22000 **Norway:** 22 67 82 53

Spain: 91 729 03 00 **Sweden:** 08 628 69 60

Switzerland: (01) 945 3242 **UK & Ireland:** (01923) 816375

31-7-95