

**Griffin & George (Sales) Limited ... : the laboratory furnishers : Ealing Road, Alperton, Wembley, Middlesex.**

**Contributors**

Griffin & George (Sales)

**Publication/Creation**

Alperton : Griffin & George, 1958.

**Persistent URL**


<https://wellcomecollection.org/works/v7ncyg5n>

**License and attribution**

Conditions of use: it is possible this item is protected by copyright and/or related rights. You are free to use this item in any way that is permitted by the copyright and related rights legislation that applies to your use. For other uses you need to obtain permission from the rights-holder(s).



Wellcome Collection  
183 Euston Road  
London NW1 2BE UK  
T +44 (0)20 7611 8722  
E [library@wellcomecollection.org](mailto:library@wellcomecollection.org)  
<https://wellcomecollection.org>

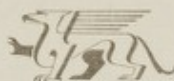

EALING ROAD, ALPERTON,  
WEMBLEY, MIDDLESEX

## GRIFFIN & GEORGE (SALES) LIMITED

Directors: K. G. SINCLAIR (Chairman), F. A. RENN (Managing), P. St. V. TABBERNER, B.A., H. C. MAYER, A.A.C.C.A. (Secretary)

THE LABORATORY FURNISHERS

TELEPHONE: PERIVALE 3344

TELEGRAMS: GRIFGEORGE, WEMBLEY

Please quote our reference:

Your reference:

FAR/P.2053

Monday, 28th July 1958.

Dear Sir,

One thousand gallons of extremely pure deionised water for little more than a shilling.

The Griffin-Raleigh Water Deioniser described in the accompanying leaflet makes use of a new process enabling mixing and reclassification of the resins to be carried out in columns down to 3 inches in diameter, thus making it a simple matter to regenerate the resins in situ. The cost of pure water is only 12s.6d. per 1,000 gallons when the apparatus is used on the relatively hard London mains supply; in areas such as Belfast and Plymouth where the raw water contains a very small proportion of dissolved electrolytes the cost is as low as 3d. per 200 gallons.

The instrument has been so designed that the methods employed for its regeneration are technically identical with those used in the regeneration of full scale industrial equipment. The diagram on the centre page of the leaflet shows the regeneration process in detail.

Two laboratory models of the Deioniser are available, carrying cylinders of 3½ and 6 inches diameter, giving 16 and 64 gallons per regeneration respectively, but we are also able to quote for larger installations as required.

We shall be very happy indeed to arrange a demonstration at your laboratory and would welcome hearing from you on this point.

Yours faithfully,  
GRIFFIN & GEORGE (SALES) LIMITED

F. A. Renn  
Managing Director

A MEMBER OF THE GRIFFIN & GEORGE GROUP OF COMPANIES  
LONDON . BIRMINGHAM . MANCHESTER . GLASGOW . EDINBURGH

	L	S
Lincoln	25	43

Lindsey	73	142
---------	----	-----

Kesteven	39	70
----------	----	----

Holland	32	61	169	316
---------	----	----	-----	-----

Retford	18	24
	<hr/> 187	<hr/> 340

M O' H	C & CB	5	
	Holland	2	
	Kesteven	4-5	
	Lindsey	10	
	Notts	2	
	Hogate	2	
	Parkborough	2	(28)
		<hr/> 205	

205

250 8 400